



perform® sterile alcohol EP

Sterile, alcohol-based ready to use solution
for disinfecting surfaces in cleanrooms



- Fast, effective and residue free
- Sterile (γ-irradiated) and microbial filtrated (0,2µm)
- Suitable for cleanrooms grade A/B, ISO 5/6
- Ready to use in a handy, sterile 500ml spray system
- Color and perfume free
- Efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms. Especially for Grade A and B clean areas, according to:

- EU Guide to Good Manufacturing Practice for medicinal products Annex 1: Manufacture of sterile medicinal products, March 2009.
- FDA Draft Guidance for Industry: Sterile drug products produced by aseptic processing – current Good Manufacturing practice; U.S. Department of Health and Human Services, Food and Drug Administration, Sept. 2004

Composition

100g contain: 25g Ethanol (94%), 35g Propan-1-ol.

Microbiological efficacy:

Efficacy:	Concentration:	Contact time:
Bactericidal, EN1276, EN13697	Ready-to-use	1 minute
Fungicidal, EN1650, EN13697		1 minute
Enveloped viruses (incl. HIV, HBV and HCV), in accordance with (German registered association for combating viral diseases)DVV-/RKI guideline		30 sec.

Chemical-physical data:

Density : appr. 0,89 g/cm³ / 20°C
Color : Colorless
pH value : appr. 7 / 20°C (undiluted)
Flash point : 27°C / Method: DIN 51755 Part 1
Form : Liquid



Special advice

Use disinfectants safely. Always read the label and product information before use.

The sterile 500ml spray system offers the following advantages:

- Convenient size for handling and storage
- Ergonomically shaped spray pump for ease of use
- Remains sterile after opening

The bottles are sterile (with irradiation indicator) and double wrapped individually. perform® sterile alcohol EP comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

Instructions for use

Apply perform® sterile alcohol EP undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. perform® sterile alcohol EP can be used for glove disinfection. Maximum application rate 50 ml/m². Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass.

Packaging unit : 10 x 500 ml bottle

Art.-no. : RREIN-PER-STEP

hans j. michael gmbh

Gewerbegebiet Hart 11, D-71554 Weissach i.T., Phone: 0049-7191-9105-0, Fax: 0049-7191-9105-19
e-mail: office@hjm-reinraum.de, website: <http://www.cleanroom-hjm.com>