



perform[®] sterile concentrate QB

Sterile, liquid concentrate



- Fast and effective
- Sterile (γ -irradiated) and microbially filtered (0,2 μ m)
- Suitable for cleanroom class A/B
- Single use bottle
- Colour and perfume free
- Efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment/devices in cleanrooms. Especially for Grade A and B clean areas, according to

- EU Guidelines to Good Manufacturing Practice for medicinal products, Annex 1: Manufacture of sterile medicinal products; March 2009
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing – current good manufacturing practice; U.S. Department of Health and Human Services, Food and Drug Administration; Sept. 2004.

Composition

100g contain: 30g Alkyl (C12-16) dimethylbenzyl ammonium chloride (ADBAC/BKC (C12-16)), 3g Polyhexamethylene biguanide (monomer: 1,5-b(trimethylen)- guanylguanidium monohydrochloride)

Microbiological efficacy:

Efficacy:	Concentration	Contact time:
Levurocidal, EN 1276, EN13697	0,5% (5ml/l)	5 min.
Enveloped viruses (incl. HIV, HBV, HCV) <i>in accordance with DVV (German registered Association for combating viral diseases) / RKI Guideline</i>	0,5% (5ml/l)	60 min.
	1% (10ml/l)	5 min.

Special advice

Use biocides safely. Always read the label and product information before use.

For all of the related safety information please refer to the Safety Data Sheet.

The practical 40ml single-use bottles are easy to open without having to use a device.

They are stable and easy to empty. Excellent material compatibility!



Instructions for use

To prepare the active solution, fill a tub or bucket with water of an appropriate quality, add the correct volume of perform sterile concentrate QB and stir gently.

For example: 1 bottle (40ml) to 8 l water = 0,5% solution

If necessary, rinse surfaces that have come in contact with the product.

Packaging unit: 20 x 40 ml

Art.-no.: RREIN-PER-STCQB

hans j. michael gmbh

Gewerbegebiet Hart 11, D-71554 Weissach i.T., Phone: 0049-7191-9105-0, Fax: 0049-7191-9105-19
e-mail: office@hjm-reinraum.de, website: <http://www.cleanroom-hjm.com>